



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


# Development of Fungal Mock Community Standards for Mycobiome Studies

Juan Lopera, PhD  
Scientist, ATCC

ASM Microbe 2019  
San Francisco, CA  
June 21, 2019

Credible Leads to Incredible™

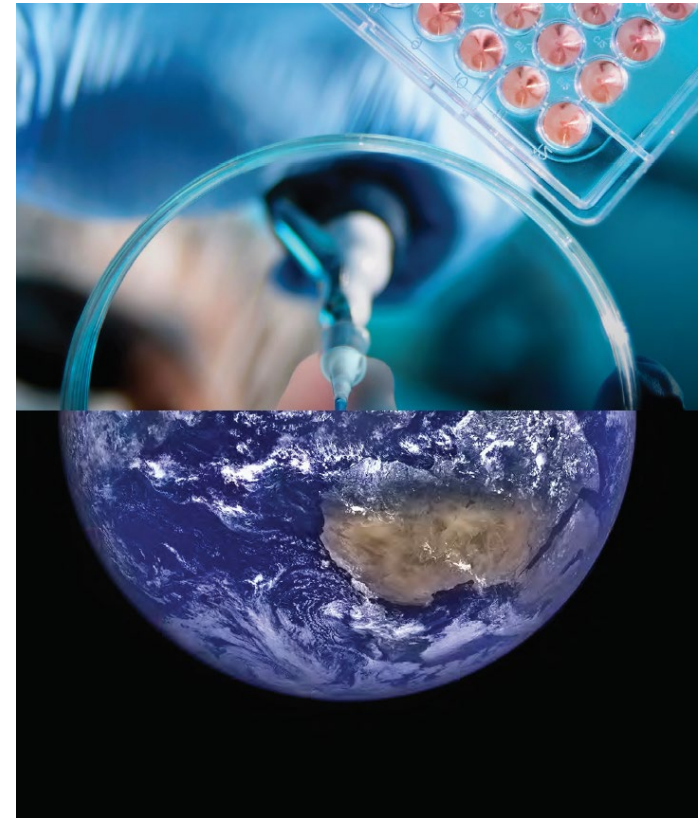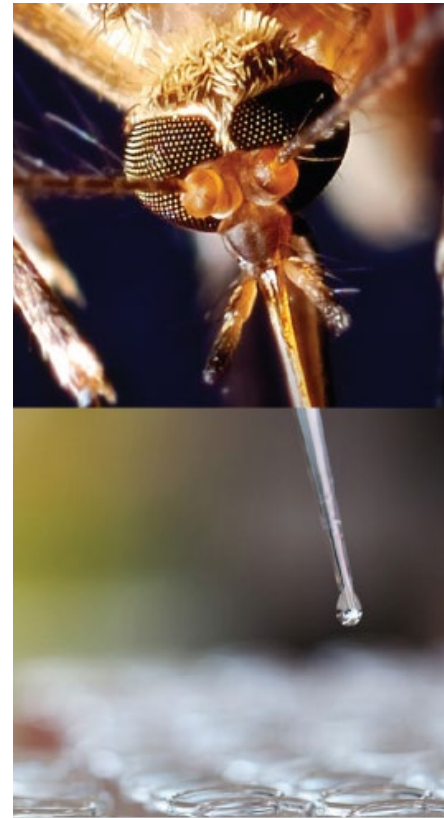

# About ATCC

*Founded in 1925, ATCC is a non-profit organization with HQ in Manassas, VA, and an R&D and Services center in Gaithersburg, MD*

*World's premier biological materials resource and standards development organization*

- 5,000 cell lines
- 80,000 microorganisms
- Genomic & synthetic nucleic acids
- Media/reagents

- ATCC collaborates with and supports the scientific community with industry-standard biological products and innovative solutions
- Growing portfolio of products and services
- Sales and distribution in 150 countries, 17 international distributors
- Talented team of 450+ employees, over one-third with advanced degrees

# An Emerging Need For Reference Materials in Mycobiome Research

frontiers in  
**MICROBIOLOGY**

**REVIEW ARTICLE**  
published: 13 February 2015  
doi: 10.3389/fmicb.2015.00089



**The lung mycobiome: an emerging field of the human respiratory microbiome**

Linh D. N. Nguyen<sup>1</sup>, Eric Viscogliosi<sup>1</sup> and Laurence Delhaes<sup>1,2\*</sup>

Cui et al. *Genome Medicine* 2013, 5:63  
<http://genomemedicine.com/content/5/7/63>



**REVIEW**

**The human mycobiome in health and disease**

Lijia Cui<sup>1</sup>, Alison Morris<sup>2</sup> and Elodie Ghedin<sup>1,3\*</sup>

**Review Article**

**The mycobiome of the human urinary tract: potential roles for fungi in urology**

A. Lenore Ackerman<sup>1</sup>, David M. Underhill<sup>2</sup>

**Cell** Article

**Temporal Stability of the Human Skin Microbiome**

Graphical Abstract

Authors  
Julia Oh, Allyson L. Byrd, Morgan Park, NISC Comparative Sequencing Program, Heidi H. Kong, Julia A. Segre




International Journal of  
*Molecular Sciences*

Review

**The Fungal Mycobiome and Its Interaction with Gut Bacteria in the Host**

Qi Hui Sam<sup>1</sup>, Matthew Wook Chang<sup>2,3</sup> and Louis Yi Ann Chai<sup>1,4,\*</sup>



CellPress

**Through the Scope Darkly:  
The Gut Mycobiome Comes into Focus**

Nydiaris Hernández-Santos<sup>1</sup> and Bruce S. Klein<sup>1,2,3,\*</sup>


Cell Host & Microbe  
**Previews**

# ATCC® Mycobiome Standards

Product description, research uses, and applications



Standard	ATCC® Catalog No.	Preparation	Description
Mycobiome	MSA-2010™	Whole Cell	Even mixture of whole cells comprising 10 fungal strains (2 x 10 <sup>6</sup> cells of each organism)
	MSA-1010™	Genomic DNA	Even mixture of genomic DNA comprising 10 fungal strains (2 x 10 <sup>6</sup> genome copies of each organism)

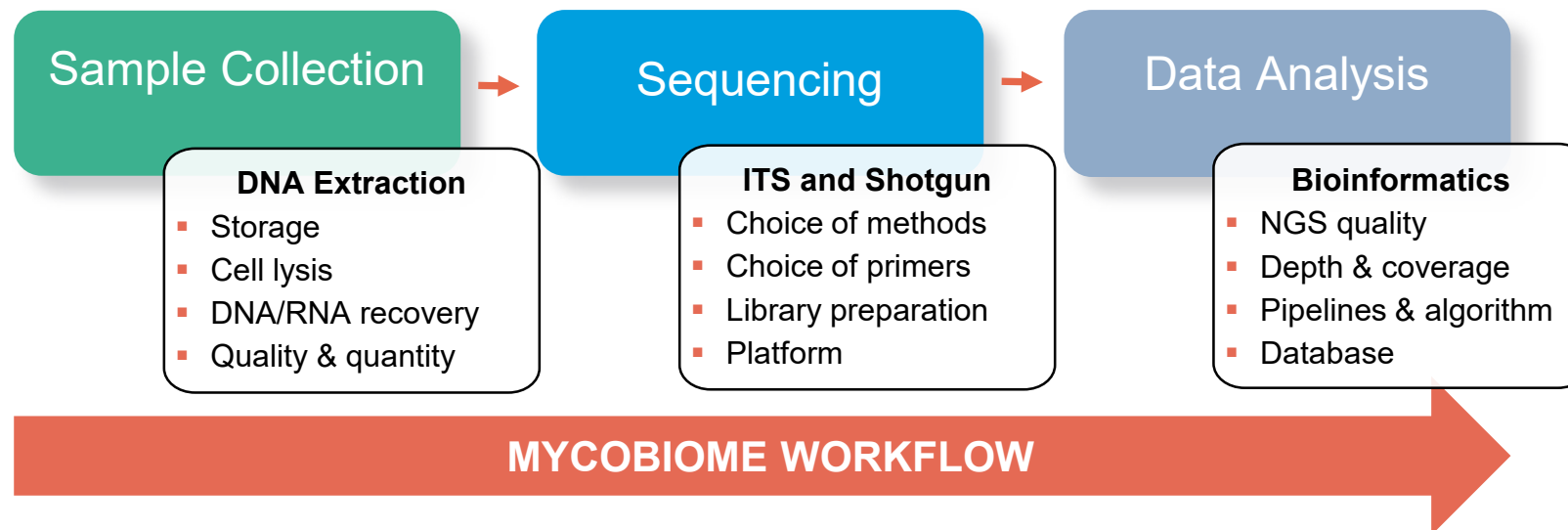

# Mycobiome Composition

Fungal strain selection, attributes, and clinical relevance

ATCC® No.	Species Name	Genome Size (Mb)	Relevancy
MYA-4609™	<i>Aspergillus fumigatus</i>	28.8	Opportunistic, airborne pathogen that is responsible for fungal infections in immunocompromised patients.
10231™	<i>Candida albicans</i>	17.1	Commensal fungus of the oral cavity that can form biofilms on denture surfaces, leading to mucosal infections.
2001™	<i>Candida glabrata</i>	12.3	Commensal fungus of the oral cavity and human gut that can acquire resistance to azole antifungals, leading to infection.
208821™	<i>Cryptococcus neoformans</i> var. <i>grubii</i>	18.9	Responsible for cryptococcal meningitis in immunosuppressed patients.
MYA-4612™	<i>Malassezia globosa</i>	9.0	Part of the normal skin flora, but can be responsible for skin diseases such as dandruff, dermatitis, and folliculitis.
201390™	<i>Saccharomyces cerevisiae</i>	12.2	Bakers' and brewers' yeast originating in food. Emerging pathogen in immunocompromised patients.
9533™	<i>Trichophyton interdigitale</i>	21.9	Can infect skin and nails to cause chronic infections such as athlete's foot and ringworm.
204094™	<i>Cutaneotrichosporon dermatis</i> ( <i>Trichosporon dermatis</i> )	23.3	Emerging opportunistic agent of invasive fungal infections, particularly in severely immunocompromised patients.
10106™	<i>Penicillium chrysogenum</i>	32.5	Spore-former, less prevalent, but can be responsible for intestinal infection in immunosuppressed patients.
36031™	<i>Fusarium keratoplasticum</i> ( <i>F. solani</i> complex)	48.6	Filamentous, opportunistic pathogen that causes fungal keratitis.


# Development of Mycobiome Standards

*Products specifications*



## Whole Cell Standards

- Authenticated ATCC cultures
- Growth and image cytometry cell counting
- Mixed in even proportion based on cell number
- Storage at 4°C

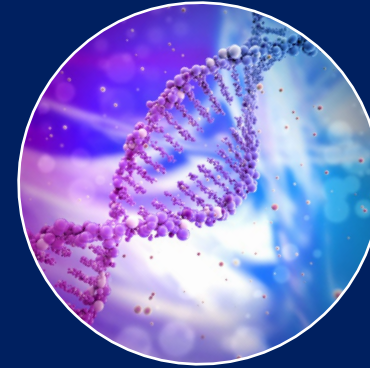

## Genomic DNA Standards


- Authenticated ATCC nucleic acids
- Fluorescent dye-based quantification
- Mixed in even proportions based on genome copy number
- Storage at -20°C

**Assay development, optimization, verification, and quality control**

# Evaluation of DNA Extraction Kits

Whole cell mock mycobiome communities can be used to evaluate the impact of DNA extraction on the mycobiome analysis


## Shotgun Metagenomic Analysis of Mycobiome Standards (ATCC® MSA-2010™ and MSA-1010™)

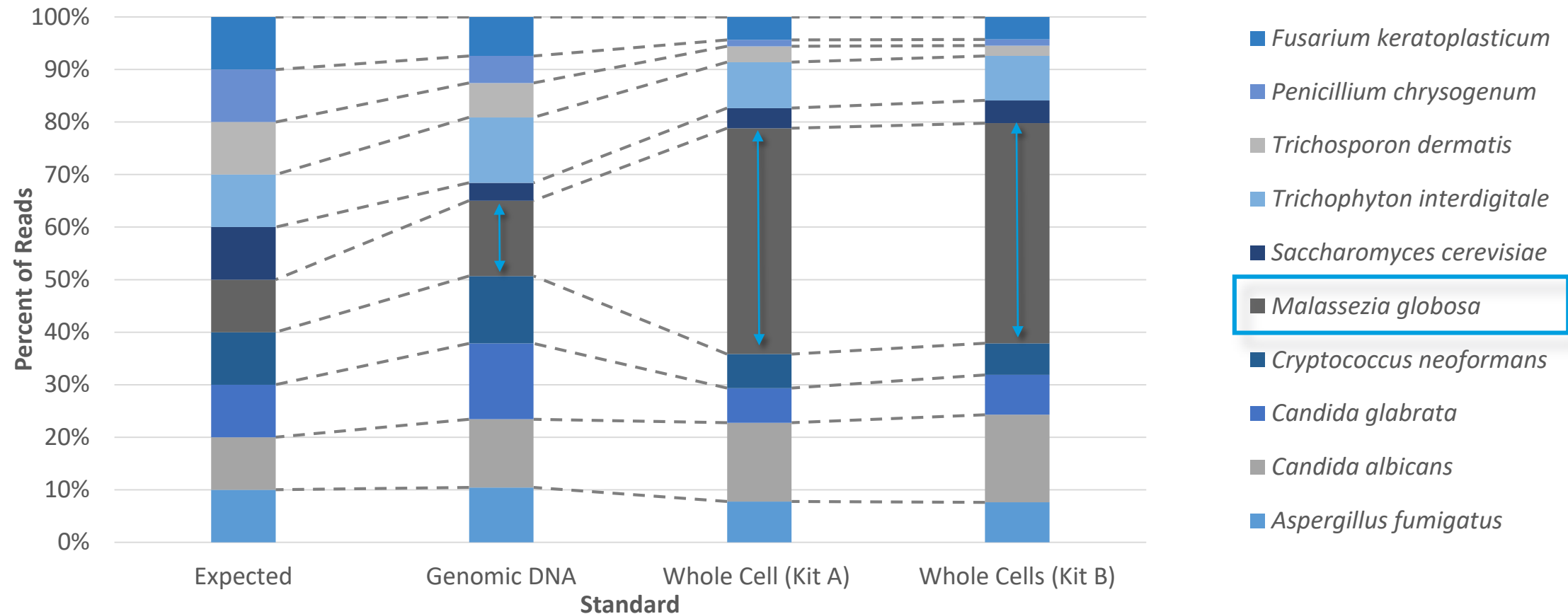


# Evaluation of ITS and Shotgun Analyses

*ITS analysis can resolve up to the genus level whereas shotgun metagenomics can identify to the species level*

**ITS Analysis – Mycobiome Standards (ATCC® MSA-1010™)**

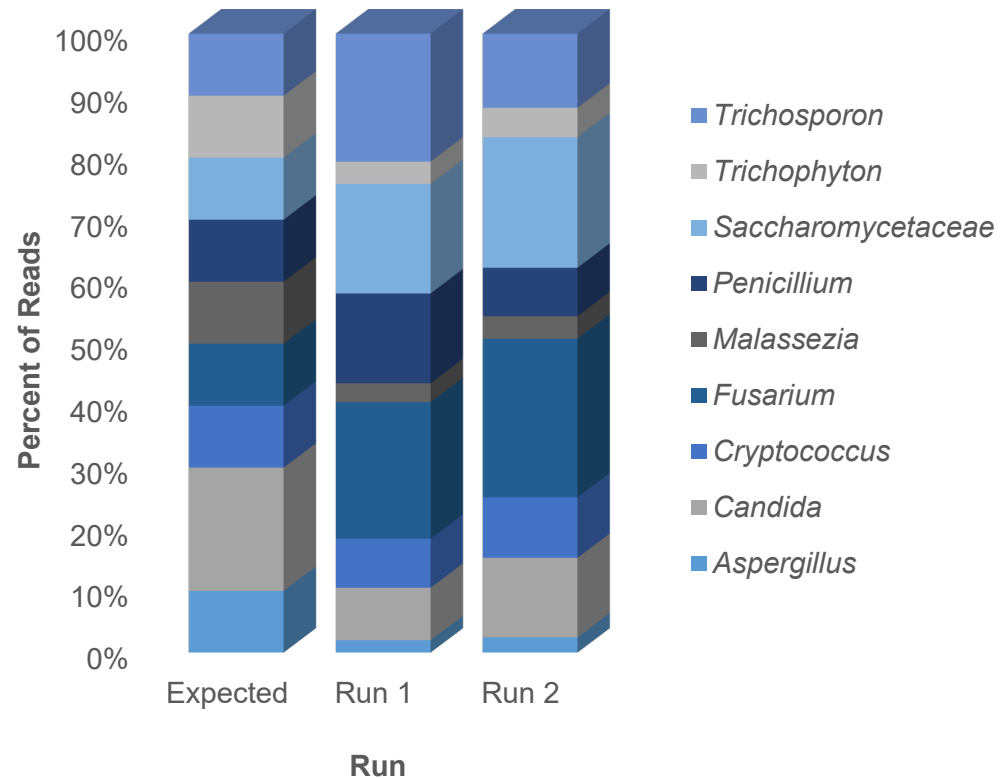

**Shotgun Analysis – Mycobiome Standards (ATCC® MSA-1010™)**

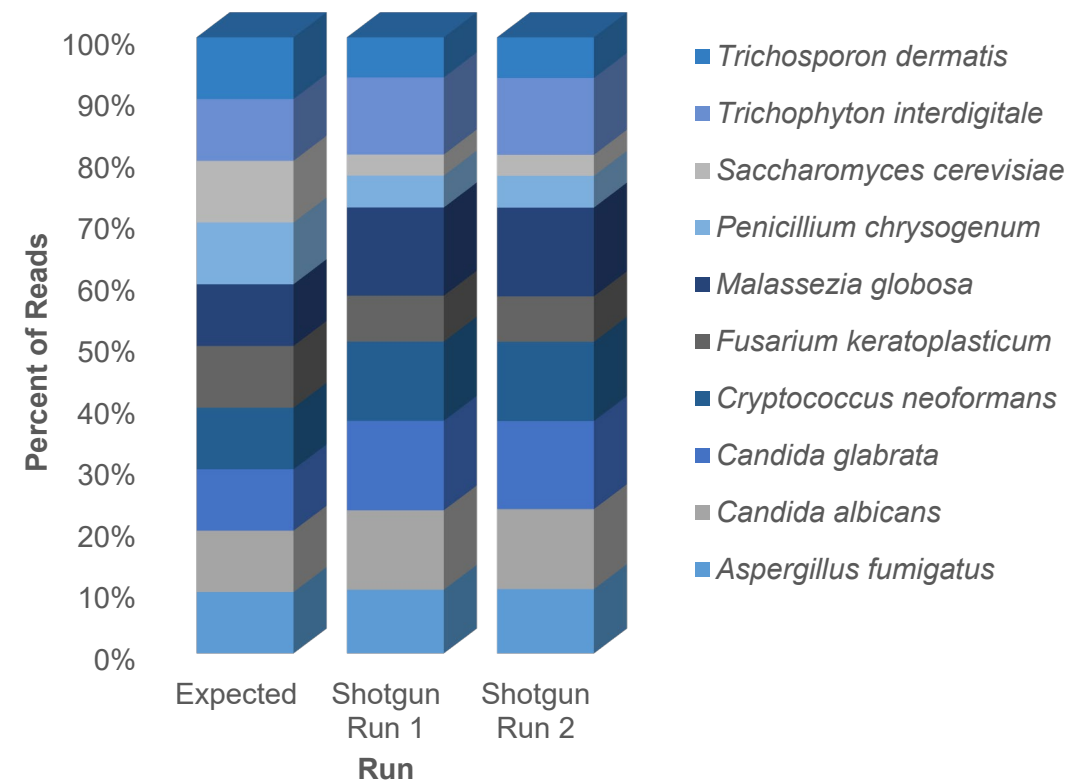




# Evaluation of Short- and Long-Read Technologies

ATCC Mycobiome Standards are platform agnostic

## Genomic DNA Standard (ATCC® MSA-1010™): Short- (Illumina) vs Long-Read (ONT) Sequencing

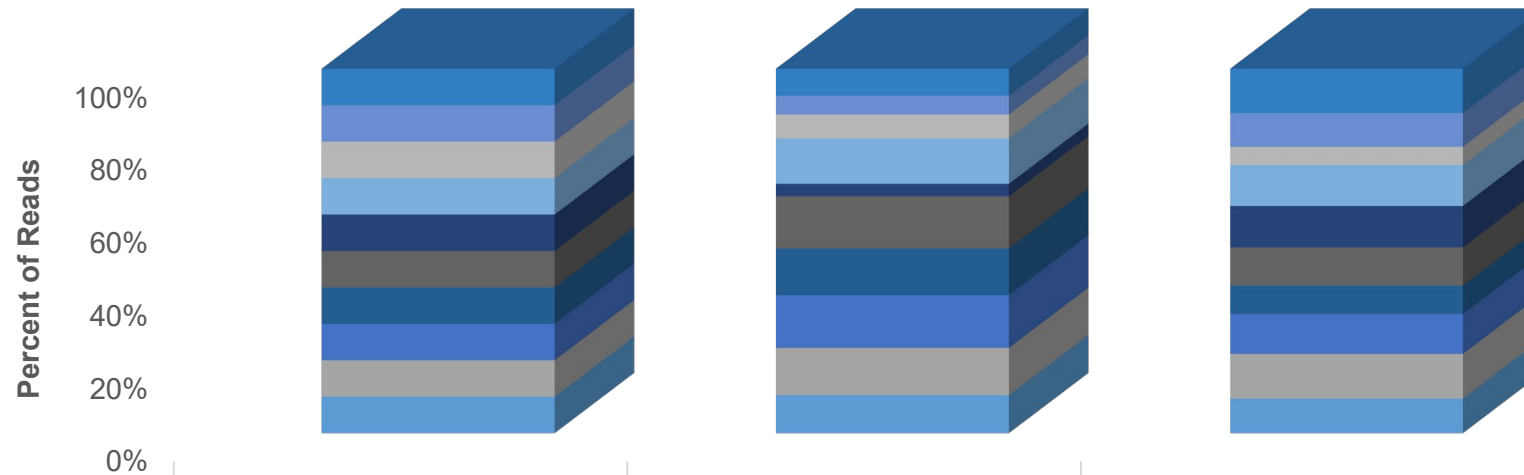


	Expected	Short-Read (Illumina)	Long-Read (ONT)
■ <i>Fusarium keratoplasticum</i>	10%	7.4%	12.2%
■ <i>Penicillium chrysogenum</i>	10%	5.2%	9.3%
■ <i>Trichosporon dermatis</i>	10%	6.5%	5.0%
■ <i>Trichophyton interdigitale</i>	10%	12.5%	11.2%
■ <i>Saccharomyces cerevisiae</i>	10%	3.4%	11.4%
■ <i>Malassezia globosa</i>	10%	14.3%	10.5%
■ <i>Cryptococcus neoformans</i>	10%	12.8%	7.7%
■ <i>Candida glabrata</i>	10%	14.4%	11.0%
■ <i>Candida albicans</i>	10%	13.0%	12.3%
■ <i>Aspergillus fumigatus</i>	10%	10.4%	9.5%

# Evaluation of Library Preparation Kits

Comparison of different library preparation kits with the genomic DNA Mycobiome Standards


## Shotgun Analysis of Genomic DNA Standards (ATCC® MSA-1010™ )

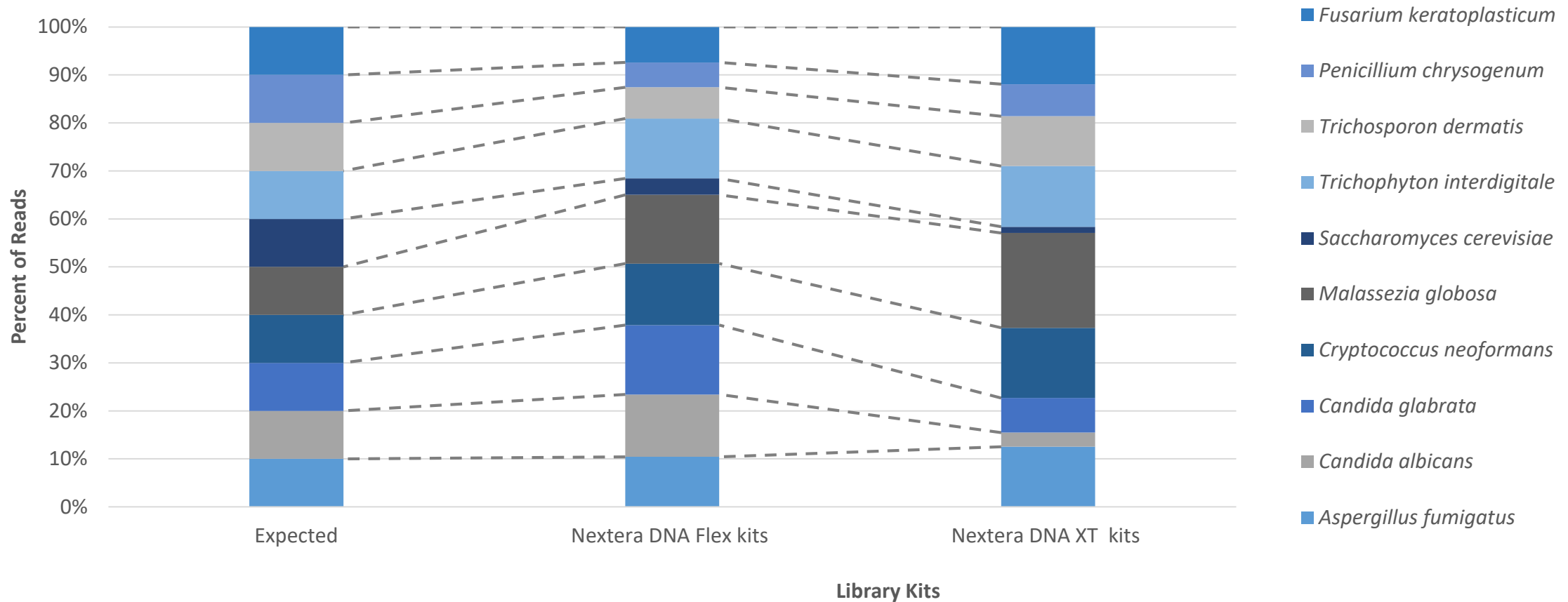

# Evaluation of Analysis Platforms and Genome Databases

Data analysis platform impacts strain identification and taxonomic resolution

## Shotgun Analysis of Genomic DNA Standards (ATCC® MSA-1010™ )

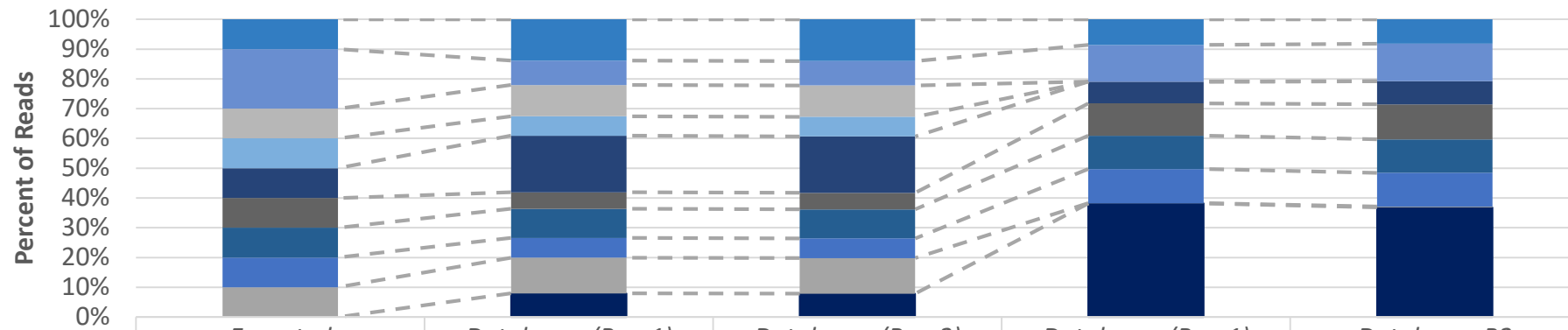

	Expected	Databases (Run 1)	Databases (Run 2)	Databases (Run 1)	Databases B2
Aspergillus	10%	13.88%	13.99%	8.60	8.16
Candida	20%	8.17%	8.22%	12.31	12.61
Cryptococcus	10%	10.52%	10.51%	0.00	0.00
Cutaneotrichosporon	10%	6.53%	6.61%	0.00	0.00
Fusarium	10%	19.00%	18.94%	7.32	7.78
Malassezia	10%	5.53%	5.55%	10.86	11.74
Penicillium	10%	9.81%	9.76%	11.17	11.25
Saccharomyces	10%	6.63%	6.62%	11.48	11.35
Trichophyton	10%	12.01%	11.98%	0.05	0.29
Unclassified	0%	7.93%	7.82%	38.21	36.82

# Acknowledgements

Monique Hunter, MS

Kendra Grosso, BS

Pushpa Gujjari, MS

Sung-Oui Suh, PhD

Anna McCluskey, BS

Stephen King, MS

Briana Benton, BS

Dev Mittar, PhD

Jonathan Hetzel, PhD, Illumina®

Nick Greenfield, MA, One Codex

© 2019 American Type Culture Collection. The ATCC trademark and trade name, and any other trademarks listed in this publication are trademarks owned by the American Type Culture Collection unless indicated otherwise. Illumina and Nextera are trademarks or registered trademarks of Illumina, Inc. Oxford Nanopore Technologies is a registered trademark of Oxford Nanopore Technologies Limited.



# Thank you!

# Questions?

Credible Leads to Incredible™

