



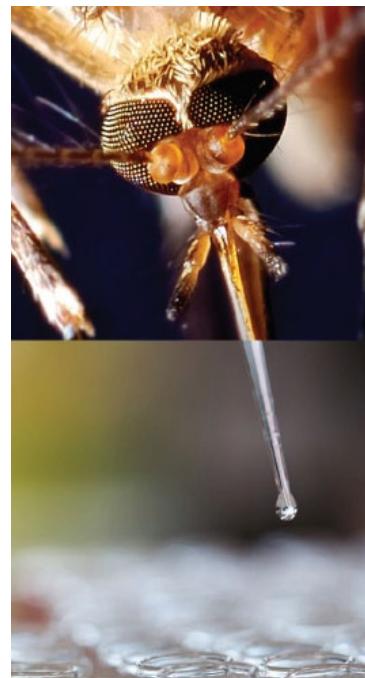

EMT Reporter Models for Cancer Research: A Window into Invasion and Metastasis

Sangeeta Kumari, MS
Senior Biologist, ATCC

Credible Leads to Incredible™

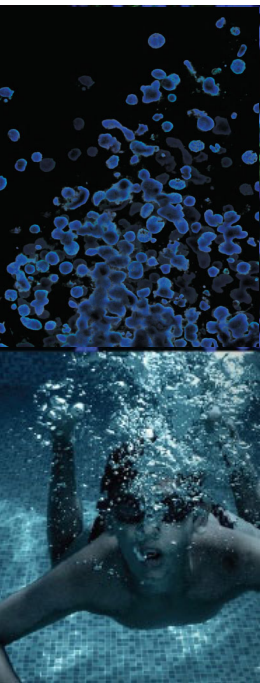


About ATCC

- Founded in 1925, ATCC is a non-profit organization with HQ in Manassas, VA, and an R&D and Services center in Gaithersburg, MD
- World's largest, most diverse biological materials and information resource for microbes – the “gold standard”
- Innovative R&D company featuring gene editing, microbiome, NGS, advanced models
- cGMP biorepository
- Partner with government, industry, and academia
- Leading global supplier of authenticated cell lines, viruses, and microbial standards
- Sales and distribution in 150 countries, 18 international distributors
- Talented team of 450+ employees, over one-third with advanced degrees

Agenda


- Background
 - EMT
 - EMT and cancer metastasis
- MCF10A ECAD EmGFP EMT
 - Gene editing strategy and confirmation
 - Morphology and growth for parental vs gene edited cells
 - Function data
- Summary

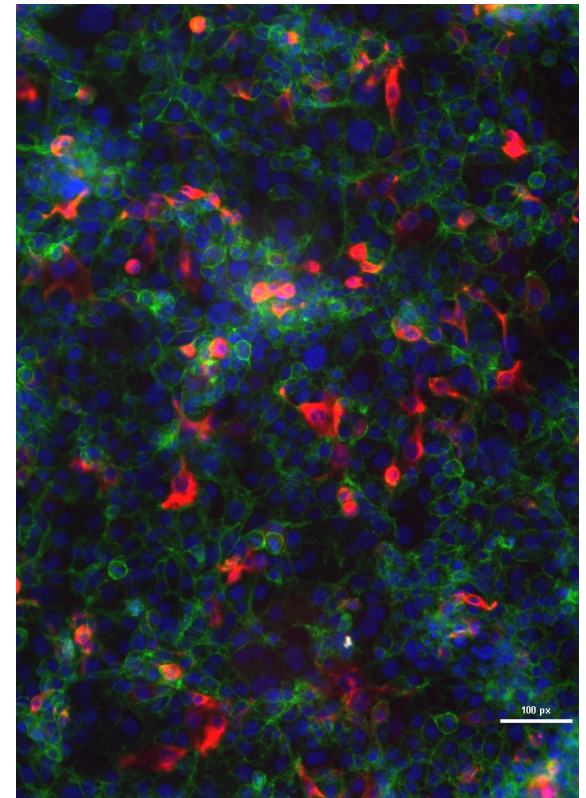

Background – EMT


- The epithelial-to-mesenchymal transition (**EMT**) is a reversible process. Epithelial cells:
 - Reduce their intercellular adhesions and proliferative capacity
 - Gain a mesenchymal phenotype with increased migratory and invasive properties

- EMT classifications and functions:
 - Implantation, embryogenesis, and organogenesis
 - Wound healing, tissue regeneration, and organ fibrosis
 - Tumor metastasis

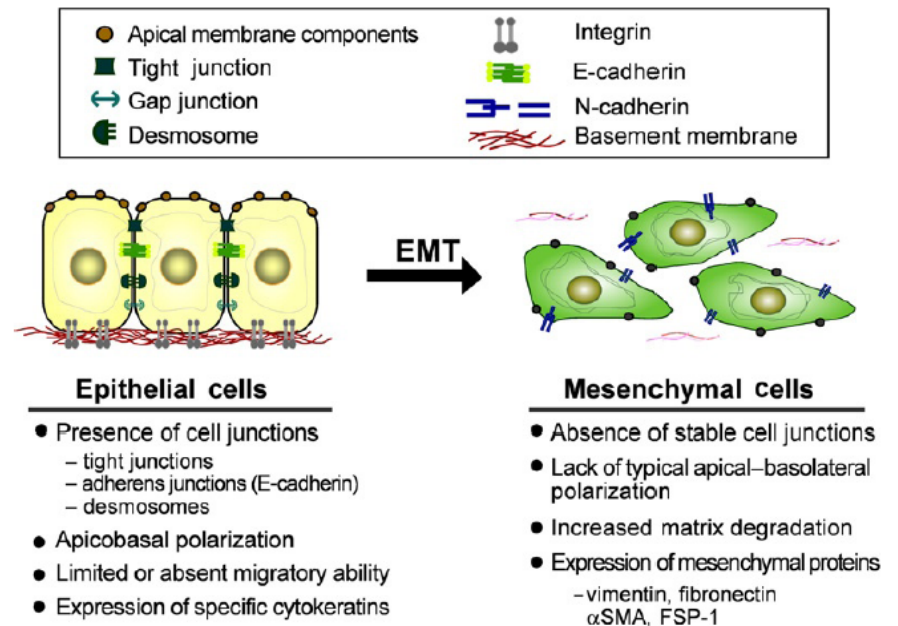

Lee et al, *International Review of Cell and Molecular Biology*, 2012.

EMT and MET in cancer progression

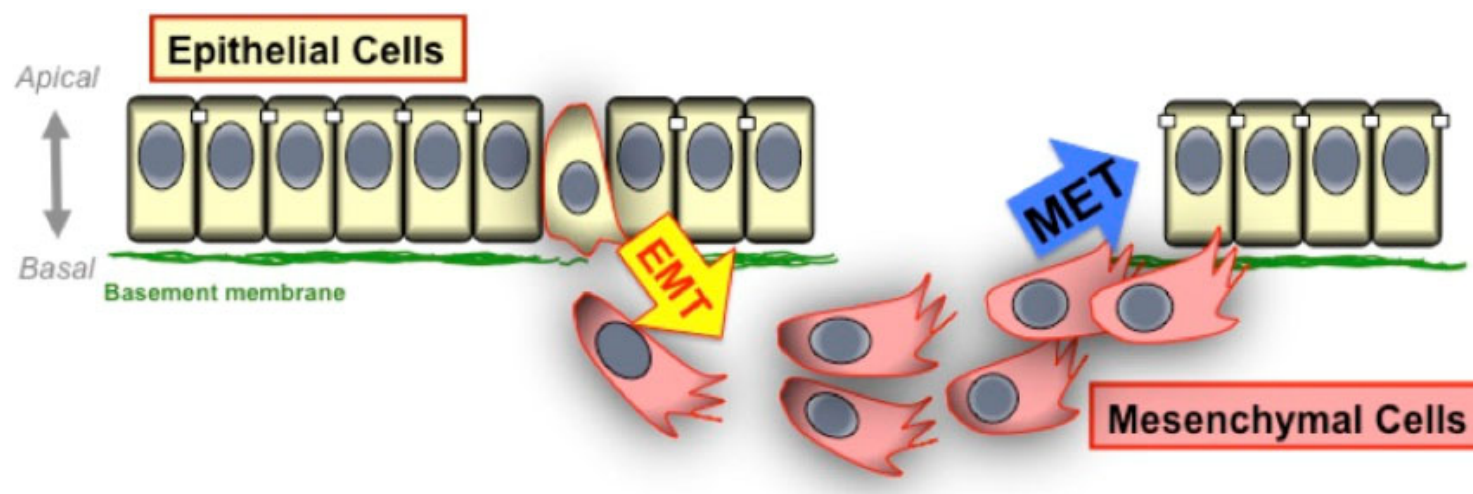


<http://murraylab.biosciences.uom.org.au/>

- The EMT process facilitates metastatic dissemination
- ATCC EMT models: **MCF10A**, **BT-474**, **A549**, **HCT116**
- ATCC MET models: **PANC-1** & **MDA-MB231**

ATCC's EMT and MET reporter cell lines

- We have developed EMT and MET reporter cell lines for use as a platform in drug screening and to learn more about the EMT/MET pathway and how it relates to cancer progression
- In these cell lines, commonly used EMT marker genes (VIM or ECAD) are tagged with a fluorescent protein to allow real-time tracking of cellular status


Designation	ATCC® No.	Tissue type/disease	EMT or MET	Marker	Availability
MCF10A ECAD EmGFP	CRL-10317EMT™	Breast epithelial cells	EMT	ECAD-GFP	Available
PANC-1 ECAD EmGFP	CRL-1649MET™	Pancreatic cancer	MET	ECAD-GFP	Available
BT-474 ECAD EmGFP	HTB-20EMT™	Breast cancer	EMT	ECAD-GFP	Available
A549 VIM RFP	CCL-185EMT™	Lung cancer	EMT	VIM-RFP	Available
HCT116 VIM RFP	CCL-247EMT™	Colorectal cancer	EMT	VIM-RFP	Available
MDA-MB-231 VIM RFP	HTB-26MET™	Breast cancer	MET	VIM-RFP	Available

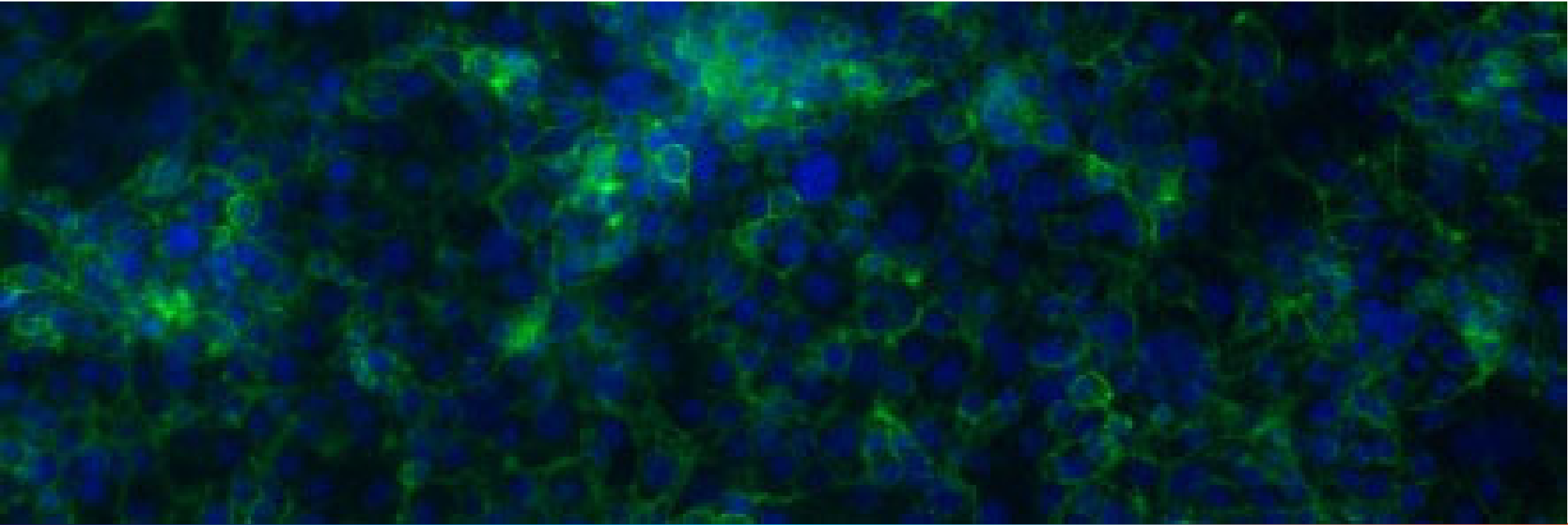

E-cadherin-EmGFP reporter lines

MCF10A ECAD EmGFP: breast epithelial cells




Generation of E-cadherin-EmGFP Knock-In Allele

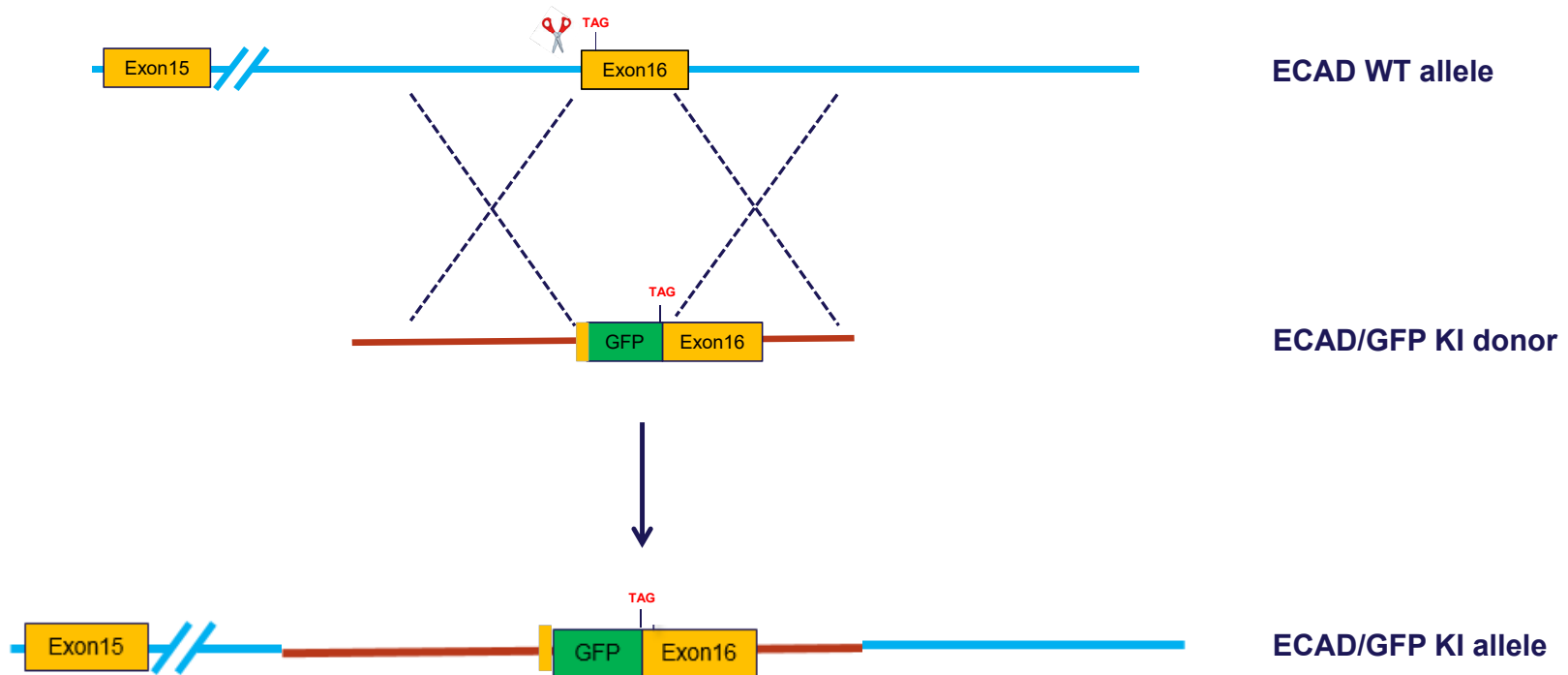


 = CRISPR/Cas9

GFP= EmGFP (Emerald Green Fluorescent Protein)

Knock-in verification at the genomic level

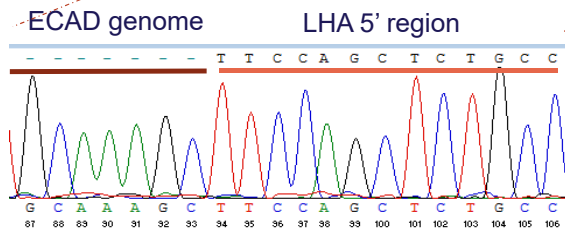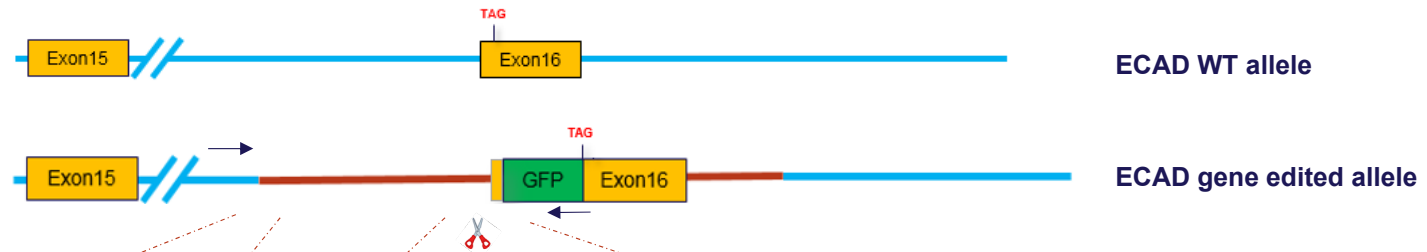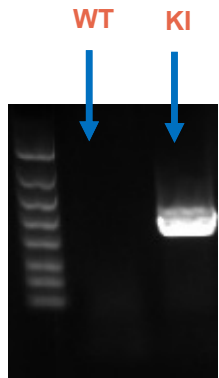


LHA junction




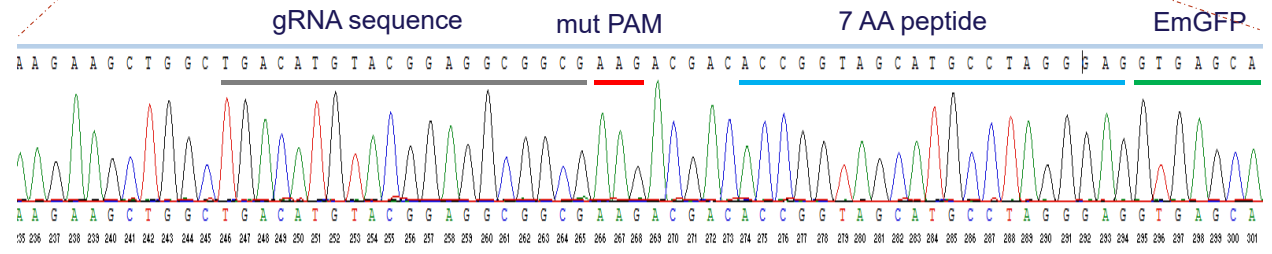


LHA/EmGFP junction

Knock-in verification at the genomic level

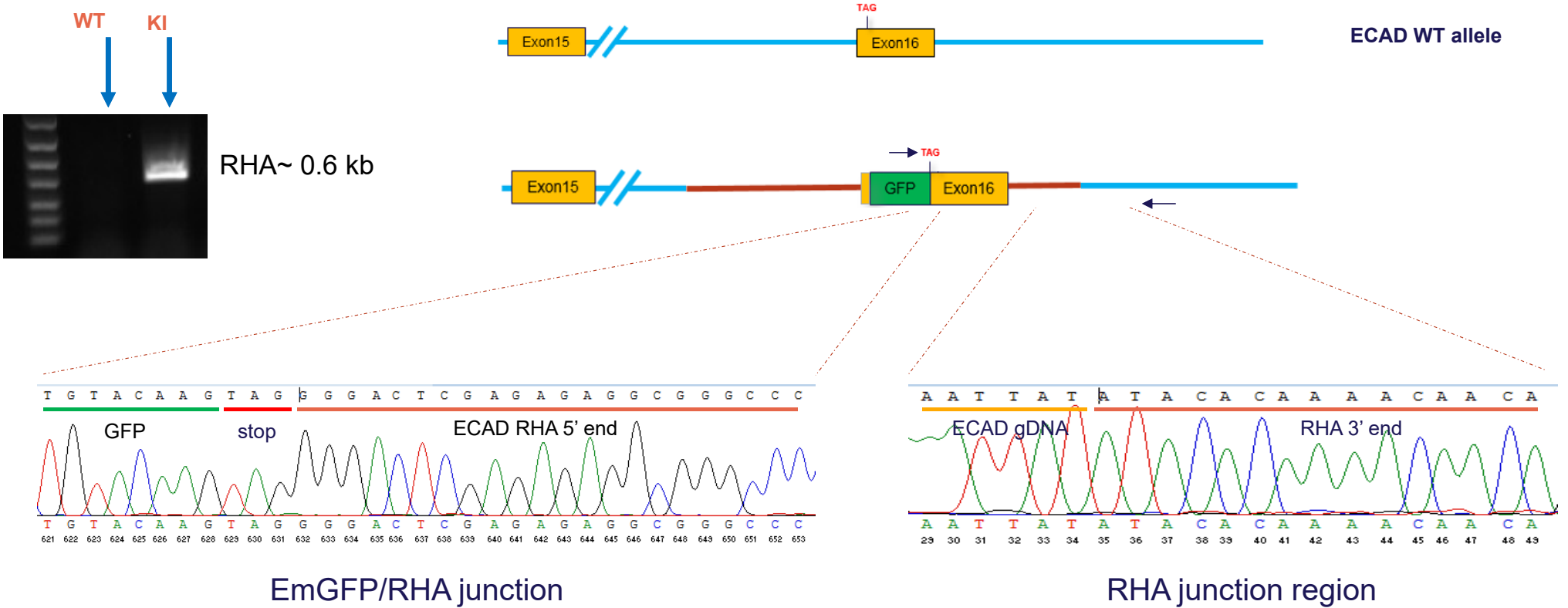

Knock-in verification at the transcriptional and translational levels

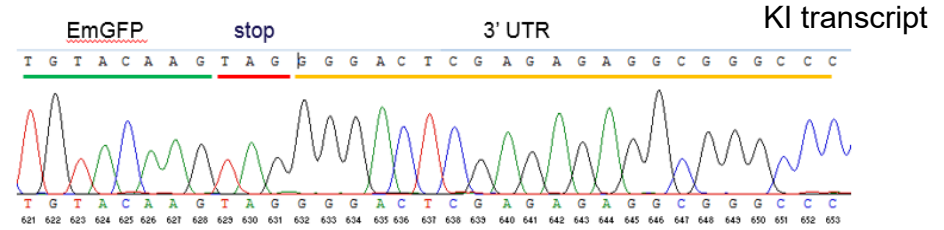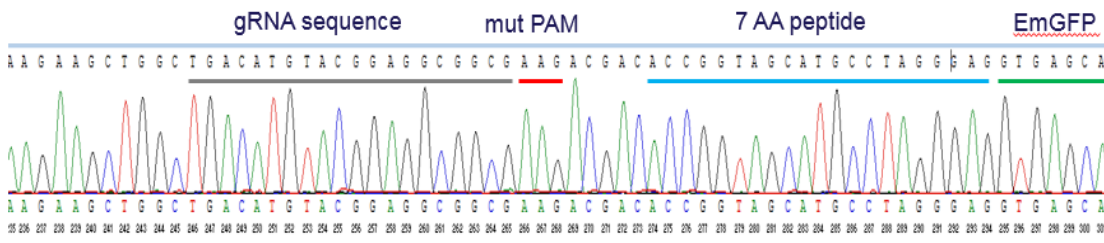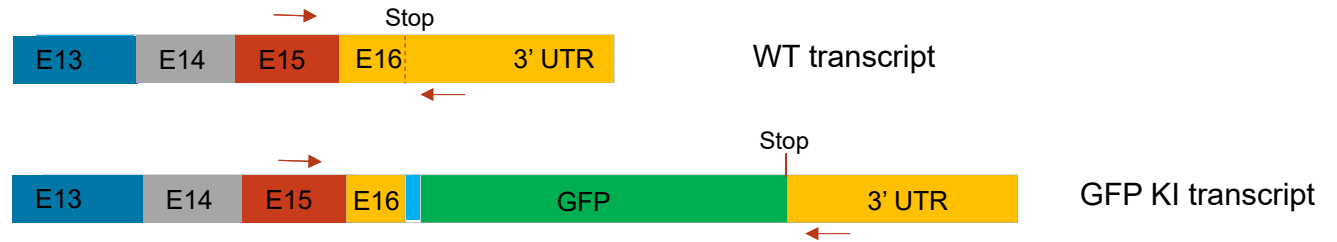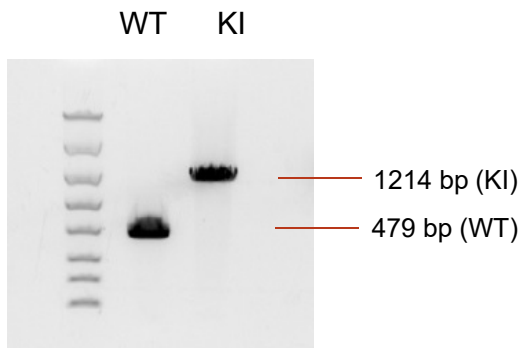


Knock-in verification at the transcriptional and translational levels

Anti-Ecadherin


Endogenous
ECAD-EmGFP

Co-localization of ECAD-GFP
expression

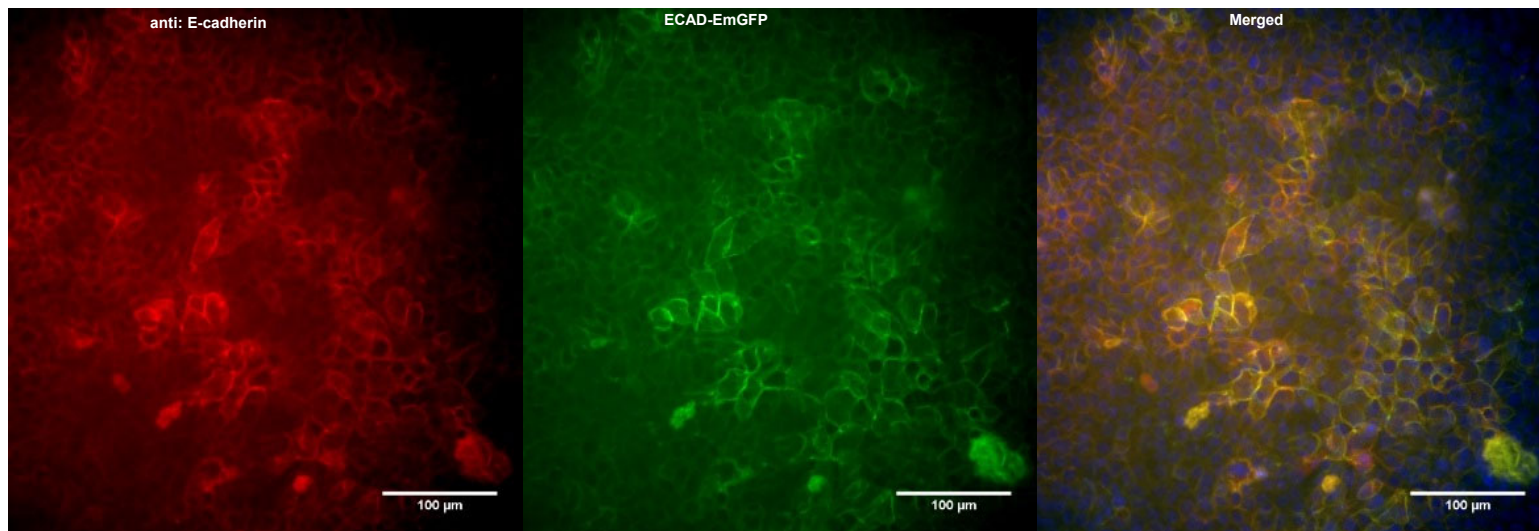


Morphology of parental and gene edited MCF 10A cells

MCF 10A WT


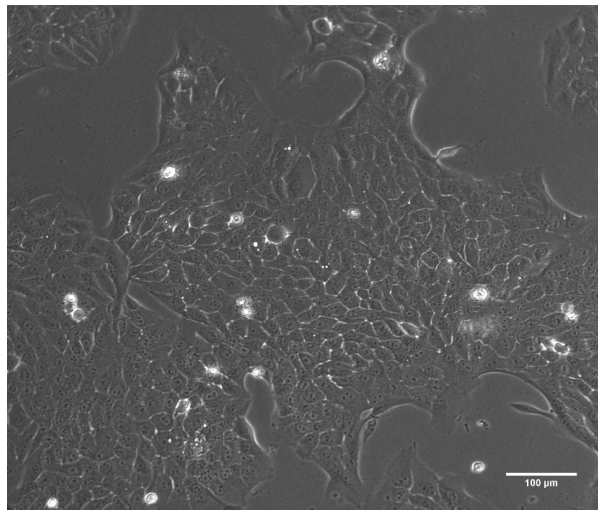


MCF 10A EMT

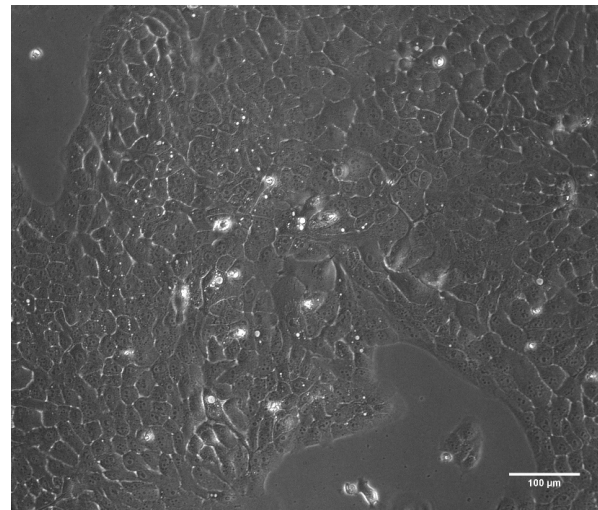

Growth media: Same as parental

Growth rate of parental and gene edited MCF 10A

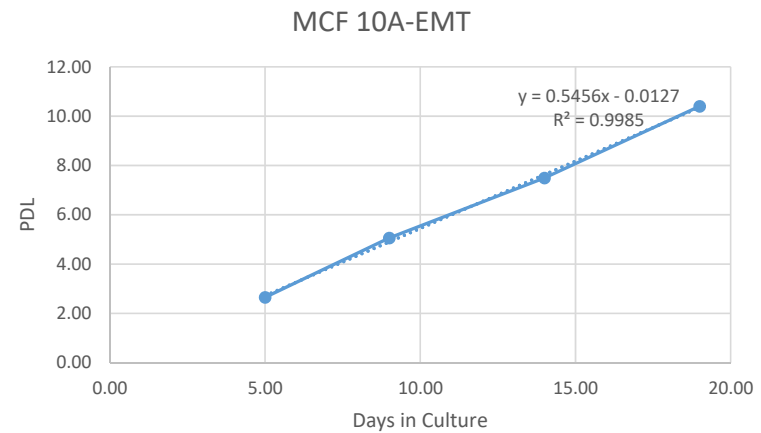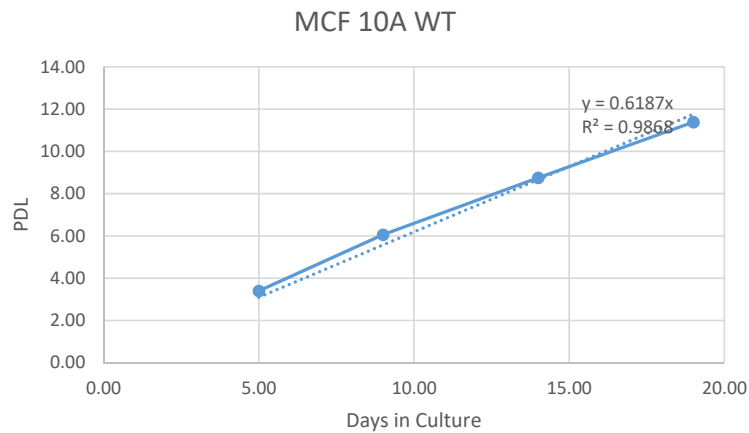


Population doubling time for MCF 10A WT = 42hrs Population doubling time for MCF 10A-EMT = 45hrs


Difference in growth = 7%

MCF 10A EMT reporter cells display morphology change upon induction


- EMT (control)

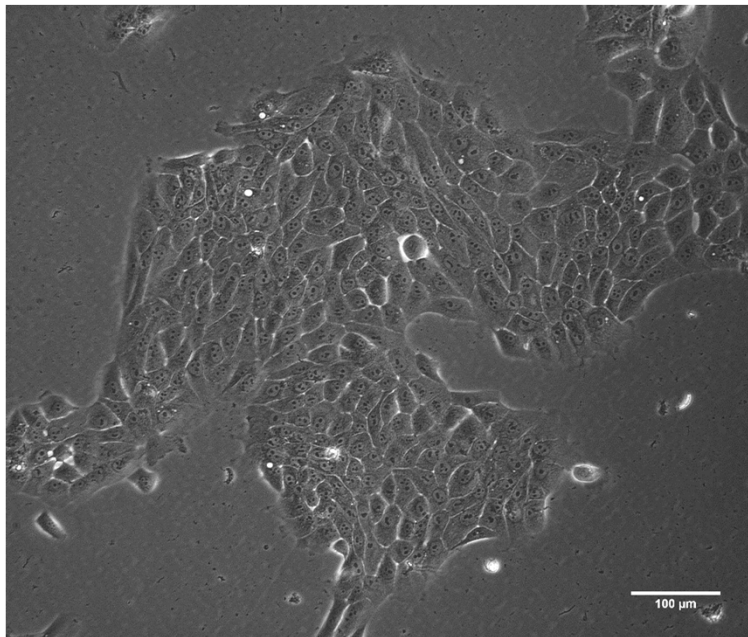



+ EMT

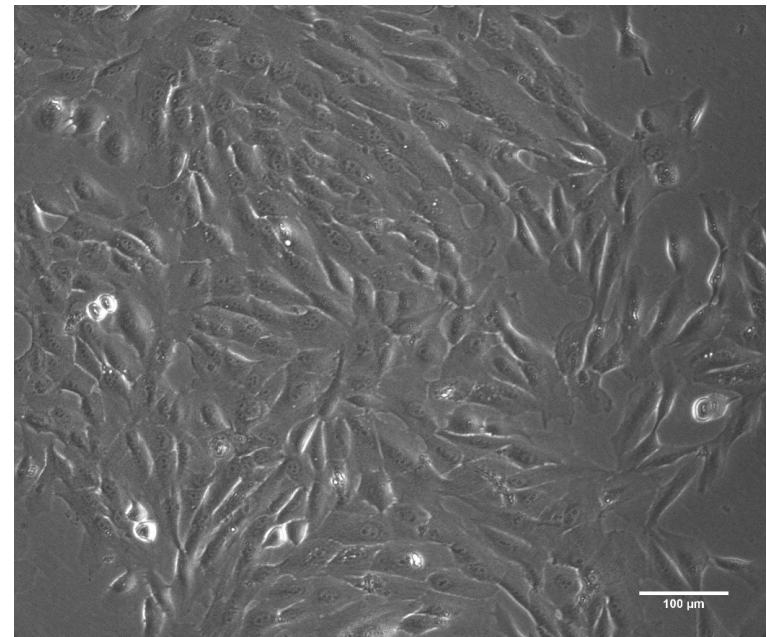



Bio-functional Data Supporting Product: Decrease in Ecad-EmGFP expression upon EMT (2 fold)

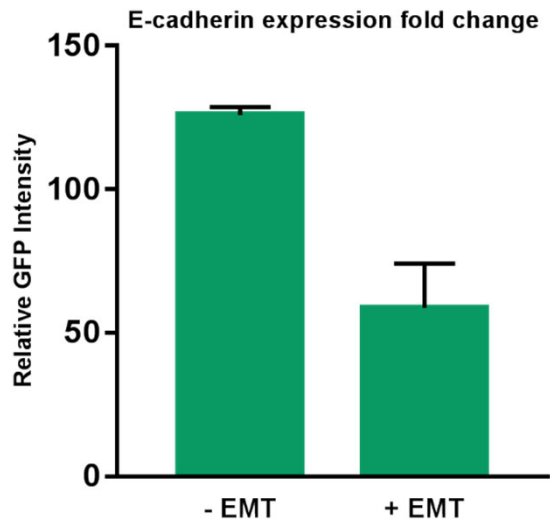




- EMT

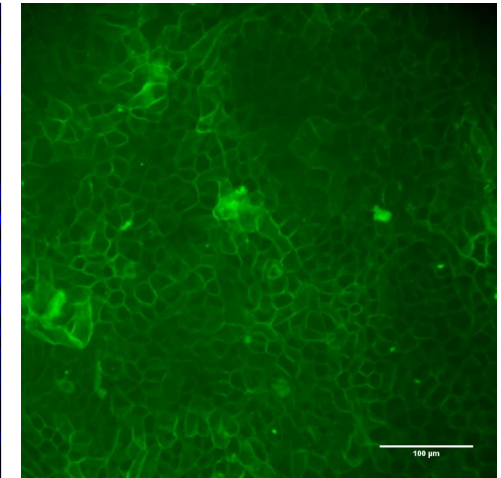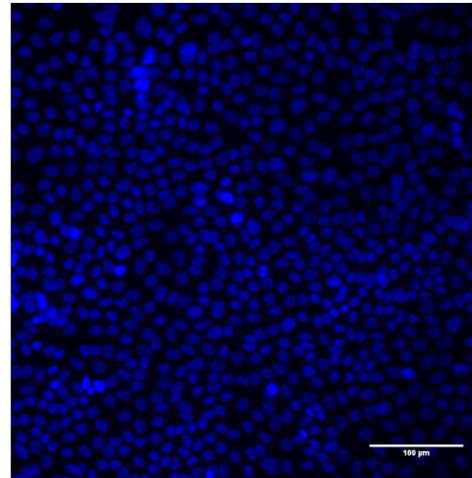


+ EMT

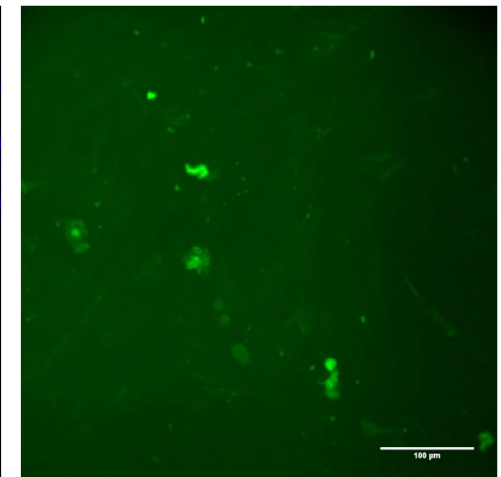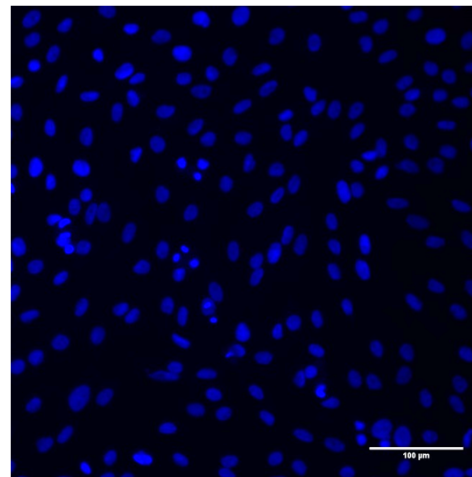


Bio-functional Data Supporting Product: ≥ 50% increase in 2nd EMT marker (Vimentin and Fibronectin) expression upon EMT

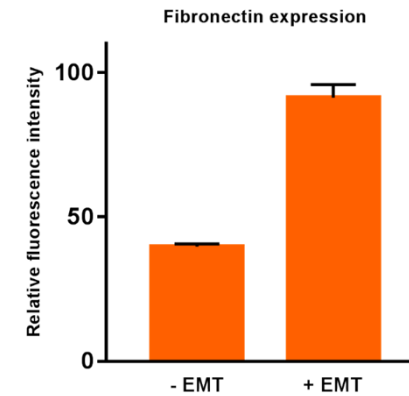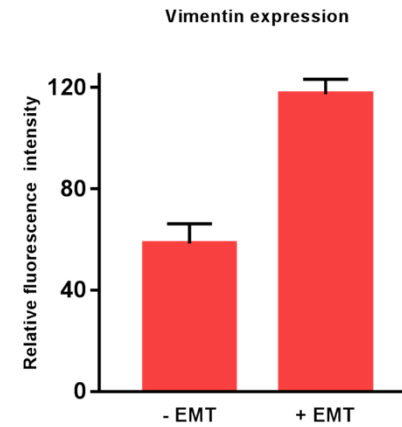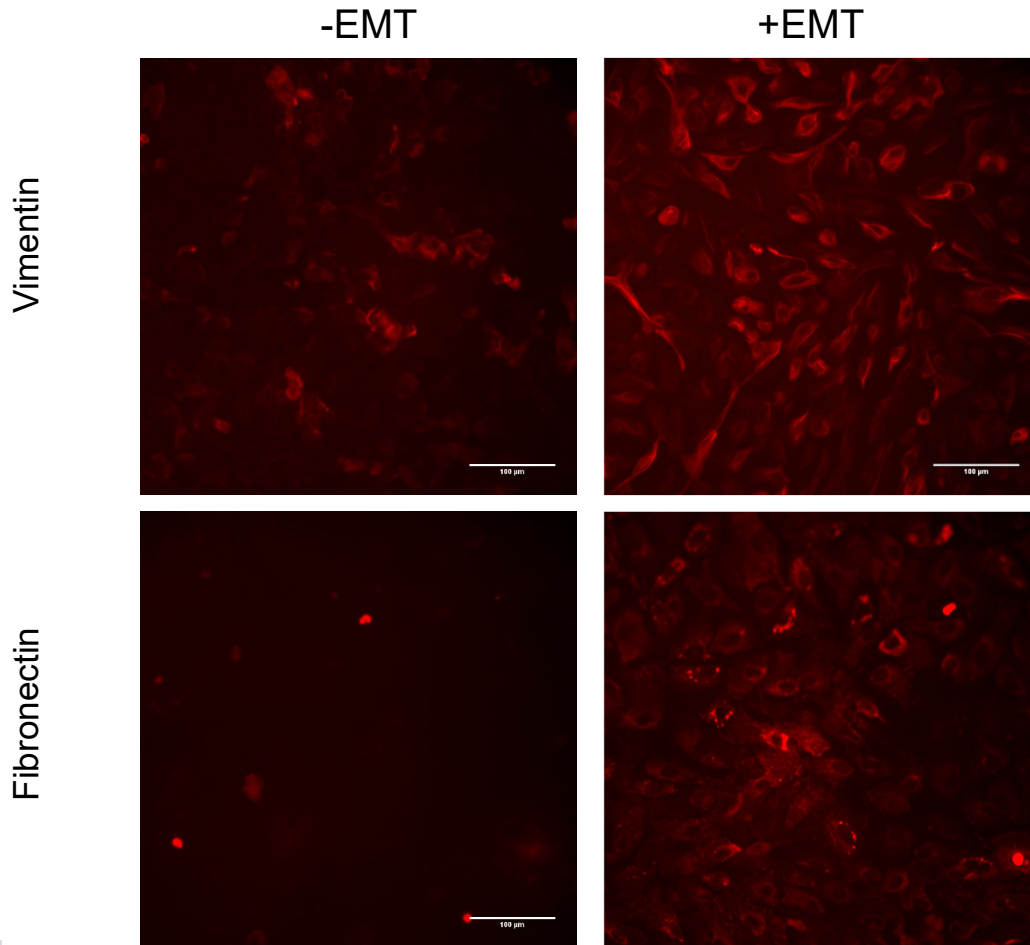


Bio-functional Data Supporting Product: MCF10A EMT cells show a significant increase in motility after EMT induction

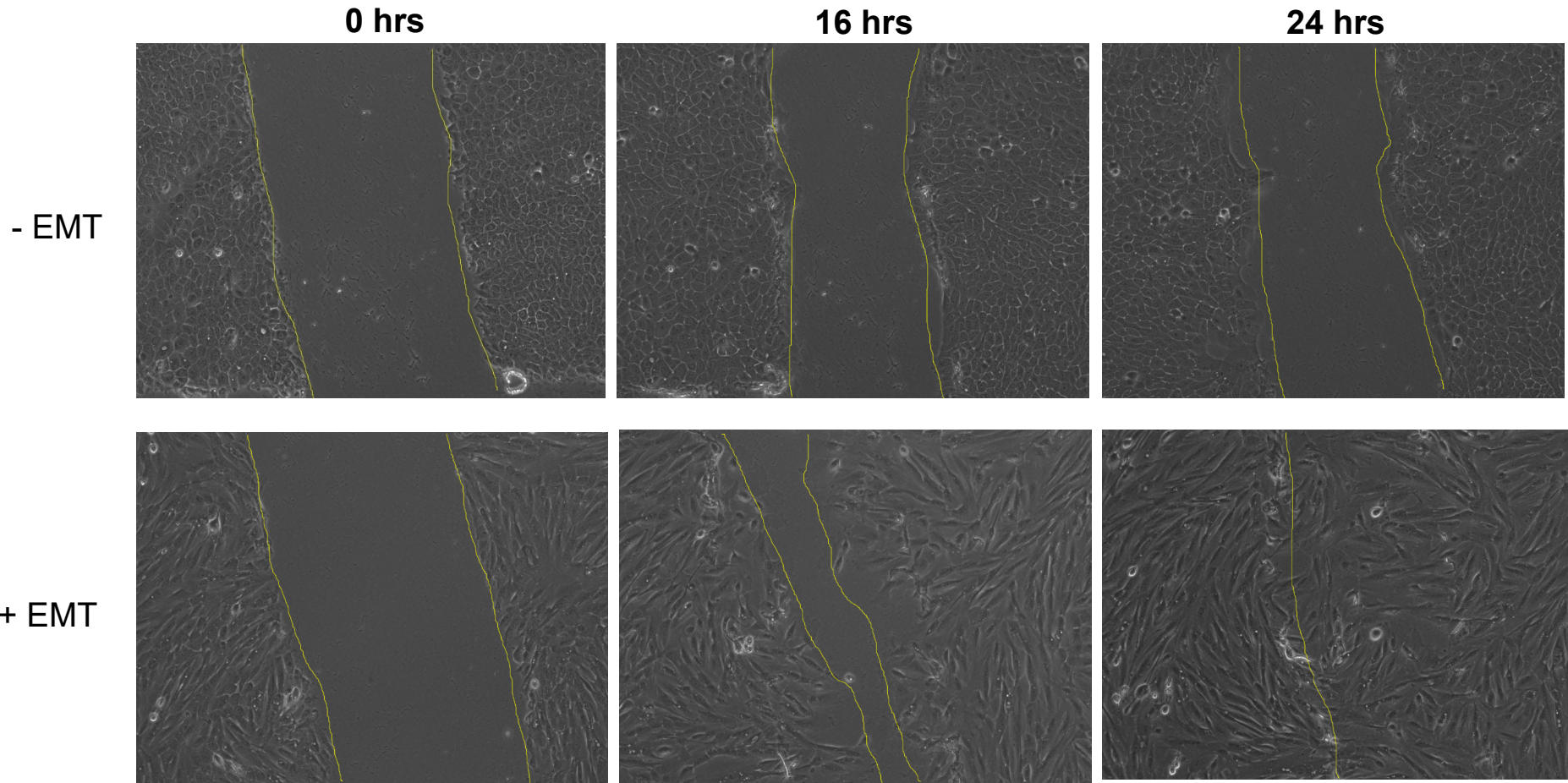

MCF10A ECAD EmGFP EMT cells show a significant increase in a migration assay

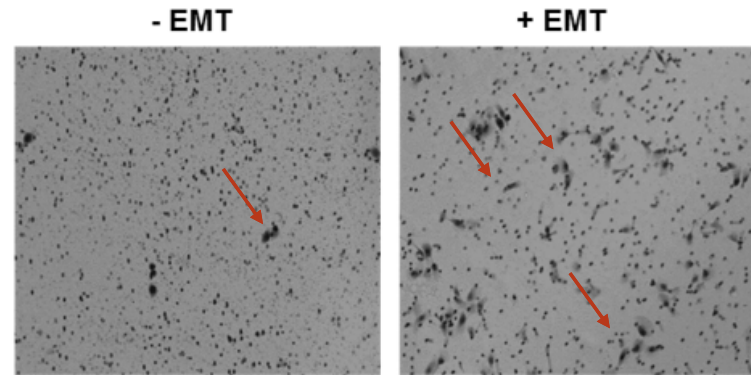


Approximately 10 fold increase in migration upon EMT induction

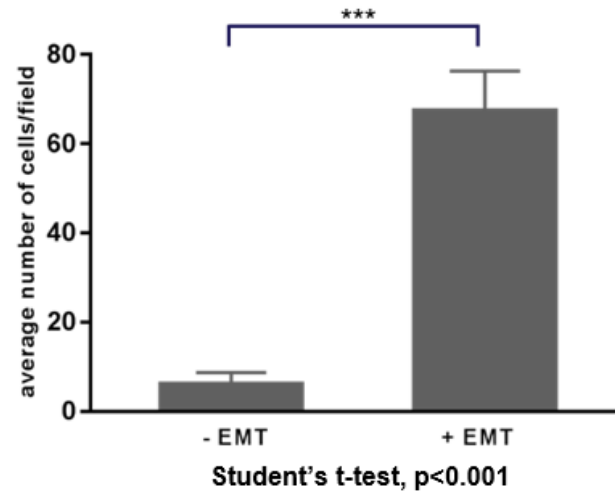

Summary

- We have created a breast epithelial, MCF 10A EMT reporter cell line
- Ecadherin, an epithelial marker is tagged to GFP using CRISPR/Cas9 gene editing technology
- Verified and validated: genomic, transcriptional, and translational levels
- Characterized with in-depth induction/transition assays
- MCF 10AEMT can be used to monitor cellular status changes in real time or as a platform to study EMT

www.atcc.org/EMT

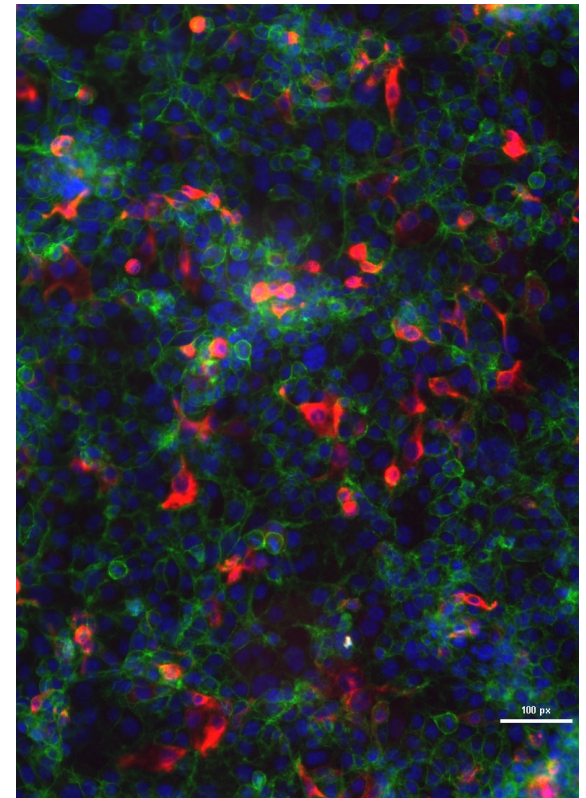

From Our Lab To Yours

Safety Tips for Returning to the Lab

To help you get back to work, we have put together a free kit containing best practices to help keep your lab infection-free:

- Poster
- Infographic
- Decals

Order the kit here:

www.atcc.org/backtothelab



© 2020 American Type Culture Collection. The ATCC trademark and trade name, and any other trademarks listed in this publication are trademarks owned by the American Type Culture Collection unless indicated otherwise.

